

# Simplifying Sample Preparation and Analysis of Released N-Glycans with a Novel Labelling Reagent Combined with HILIC Chromatography

Jennifer Fournier, Matthew A. Lauber, Sean M. McCarthy, and Stephan M. Koza  
Waters Corporation, 34 Maple Street, Milford, MA 01757 USA

The post-translational modification of proteins with oligosaccharides, known as glycosylation, can affect both the structure and function of a protein [1-3]. Not surprisingly, the glycan profile of a biopharmaceutical is commonly defined as a critical quality attribute [4-5], since it can be a measure of efficacy and immunogenicity, as well as an indicator of manufacturing conditions [6-7]. N-glycosylation of proteins, such as monoclonal antibodies (mAbs), is consequently monitored during the development and lot release of biotherapeutics. Very often these assays are performed on N-glycans after they are released from their counterpart proteins [8]. Since glycans are highly polar molecular structures, they are amenable to hydrophilic interaction liquid chromatography (HILIC) based techniques for purification and separation (Figure 1). Rudd and co-workers have clearly shown the utility of HILIC for mapping the N-glycan heterogeneity of protein samples [9-15]. In such a method, the released N-glycans may be derivatised so that they bear a chromophore or fluorophore and can be detected optically following the chromatographic separation. Unfortunately, conventional approaches to the preparation of N-glycans for HILIC analysis are either laborious or require compromises in MS or optical sensitivity [16-17]. This limits the throughput or detail by which glycosylation can be characterised or monitored. In addition, many of these methods, due to their complexity, are challenging to transfer throughout an organisation. Thus, it would be advantageous for an N-glycan sample preparation protocol to be streamlined and robust, while facilitating sensitive MS and optical detection.

## Accelerating Derivatisation

N-Glycan analysis methods often rely on an analyte derivatisation procedure that employs reductive amination. Reductive amination labels the reducing, aldehyde termini that form on N-glycans only after they hydrolyse from their glycosylamine form. This reaction requires an anhydrous reaction, which begins with a dry down step followed by a 2-3 hour chemical conversion process (Figure 2) [16-17].

These methods often contain numerous steps that require significant levels of

expertise to optimise and because of the complex and stringent nature of this chemistry there is a high potential for analyst error which can result in repeating the sample preparation. To address this challenge, rapid tagging reagents have recently been introduced that can be used in place of conventional reductive amination labels, such as 2-aminobenzamide (2AB) [18]. These reagents target the native glycosylamines which are initially formed during enzymatic release from protein, to yield highly stable urea-linked derivatives. Although these reagents have accelerated labelling and quickly introduce a fluorophore for chromatographic detection, they have lacked the chemical properties needed to facilitate mass spectral identification of low abundance N-glycans. An improvement upon this approach for labelling

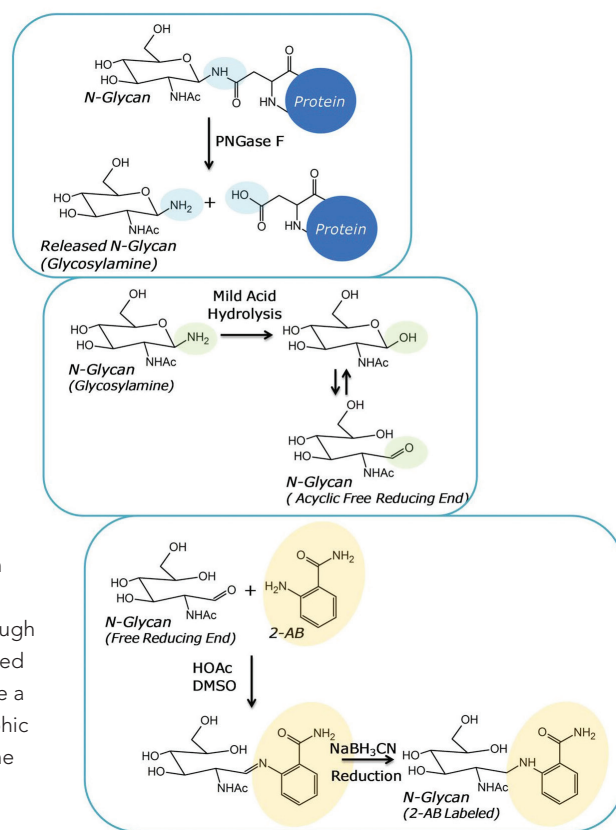


Figure 2. Reductive Amination of N-Glycans

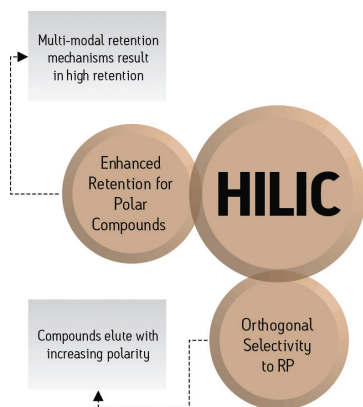


Figure 1. Benefits of HILIC SPE for Glycans.

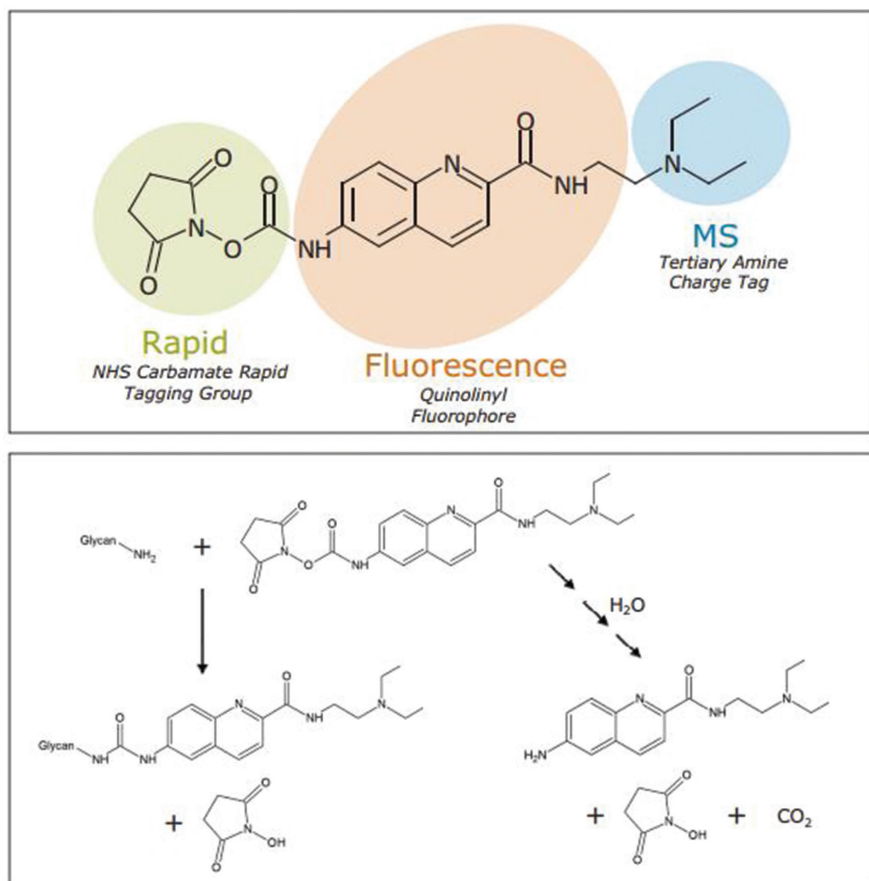


Figure 3. The chemical structure of RapiFluor-MS and its reaction with an N-glycosylamine.

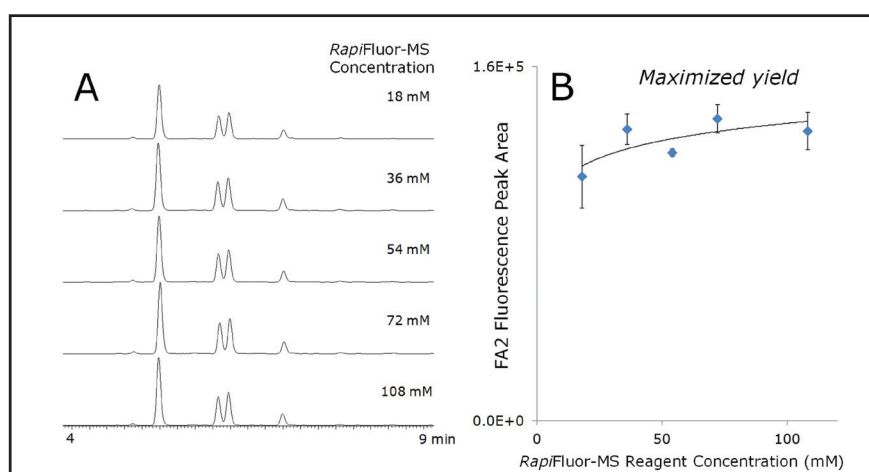


Figure 4. Optimisation of labelling reagent molar excess for RapiFluor-MS labelling (A) Fluorescence chromatograms for labelled glycans obtained by titration of 0.36 mg/mL deglycosylated mAb with varying concentrations of RapiFluor-MS reagent. Separations were performed with labelled glycans from 0.4  $\mu$ g of glycoprotein and a 2.1 x 50 mm ACQUITY UPLC BEH Amide 130Å 1.7  $\mu$ m column. (B) Fluorescence peak area as a function of RapiFluor-MS reagent concentration.

released N-glycans would be designed to also derivatise the glycosylamine with a label that would allow for enhanced fluorescence and MS detectability. **Figure 3** shows a purposefully designed compound called RapiFluor-MS™ that has these features.

In a 5 minute reaction, N-glycosylamines are labelled with RapiFluor-MS, a reagent comprised of an N-hydroxysuccinimide (NHS) carbamate rapid tagging group, a high quantum yield quinolone fluorophore,

and a basic tertiary amine that enhances ionisation efficiency. This approach is dependent on a quick deglycosylation reaction that limits glycosylamine hydrolysis, a reaction that has been determined to have a half-life of approximately 2 hours at 50°C in the pH 7.9 digest buffer used. To this end, a rapid, PNGase F deglycosylation produces complete deglycosylation of most glycoproteins in 5 minutes. Coupled with nearly quantitative labelling it is

predicted that less than 3% of the released glycosylamine will be unlabelled under the proposed conditions [19].

### Building Robustness into N-Glycan Sample Preparations

Gains in efficiency and sensitivity are important, but so is having a protocol that is robust and able to produce consistent results. Every step in a protocol can potentially introduce bias and sample loss, for example a step to deplete the protein from the sample after deglycosylation. Not performing a protein removal step results in higher and more consistent glycan recovery when a molar excess of the labelling reagent based on glycoprotein concentration is employed. Studies have been performed with RapiFluor-MS labelling to show that the use of both higher and lower than recommended reagent produce comparable fluorescent profiles, indicating a high degree of robustness in the labelling (**Figure 4**).

While a clean-up step prior to labelling is avoided in this N-glycan sample preparation method, it is advantageous to have a clean-up step for the labelled glycans prior to HILIC chromatographic analysis, so that baseline interferences are minimised, resulting in improved relative quantitation from fluorescence chromatograms. However, extraction of the labelled glycans from reaction by-products is often a step that introduces bias. As mentioned previously, HILIC is an ideal mode for retaining glycans due to their highly polar nature. Therefore, a procedure has been optimised using a HILIC SPE process, involving a silica based aminopropyl sorbent. This particular stationary phase retains polar compounds with a surface that is both strongly hydrophilic and weakly basic, which allows an analyst to take advantage of electrostatic repulsion and ion exchange to improve yield and purity of the labelled N-glycans. The charge of the stationary phase surface is dependent on pH, at higher pH values the ionisation of basic analytes and the aminopropyl ligand is reduced, however the ionisation of the surface silanols is increased. Therefore, it is critical to choose an eluent that will have a fixed stable pH each time a sample preparation is performed (**Figure 5**).

### Combining HILIC Chromatography with Mass Detection

Finally, as with some traditional labelling chemistries such as 2-AB and 2-AA,

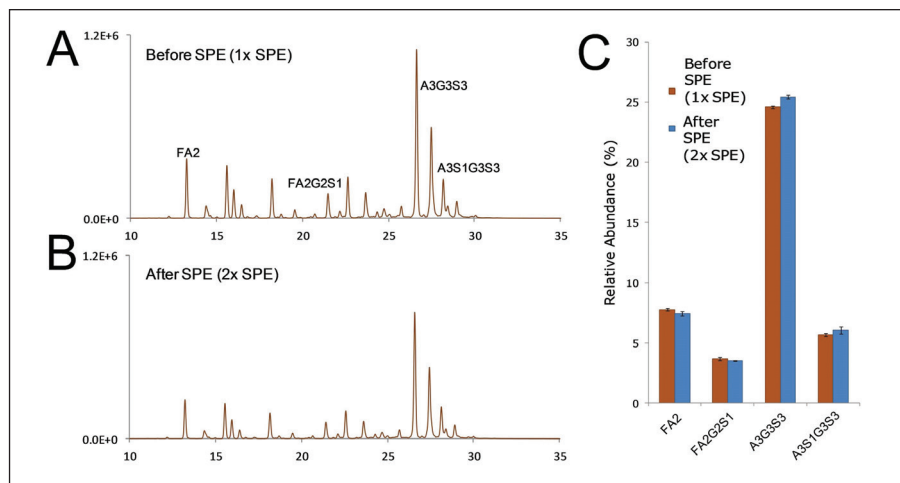


Figure 5. Extraction of RapiFluor-MS labelled N-glycans by SPE with a GlycoWorks HILIC  $\mu$ Elution plate. (A) A test mixture comprised of RapiFluor-MS labelled glycans from pooled human IgG and bovine fetuin separated on a 2.1 x 150 mm ACQUITY UPLC BEH Amide 130Å 1.7 $\mu$ m column and detected via fluorescence (Labelled N-glycans from 0.4  $\mu$ g glycoprotein, 10  $\mu$ L injection of ACN/DMF diluted sample). (B) The test mixture after extraction by HILIC SPE. (C) Relative abundances determined for a set of RapiFluor-MS labelled glycans before and after GlycoWorks HILIC SPE.

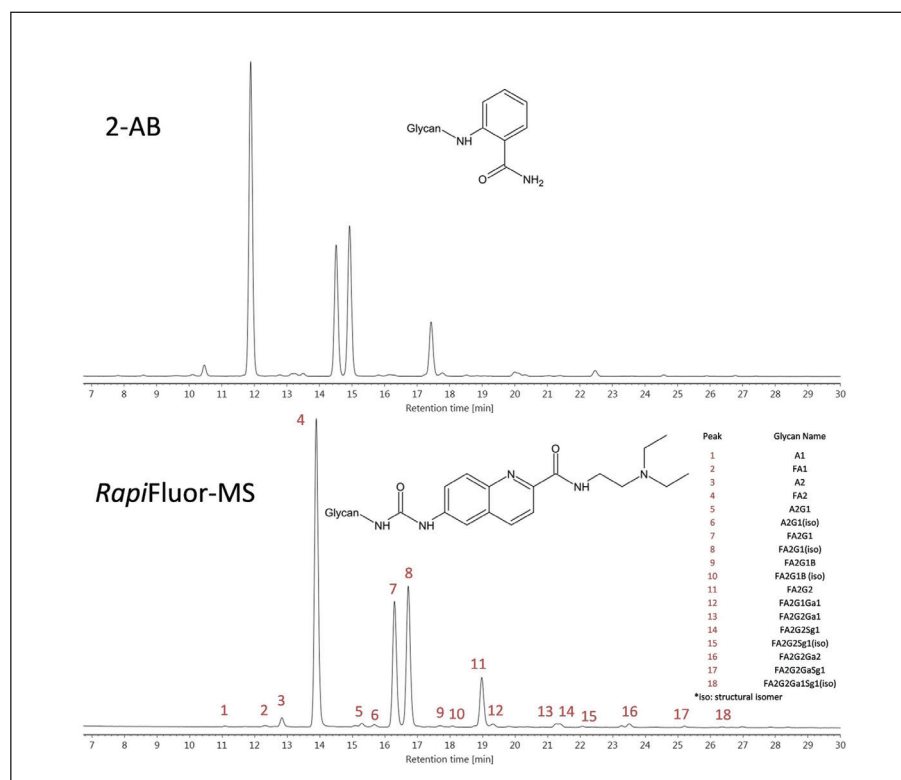


Figure 6. Similarity between 2-AB and RapiFluor-MS N-glycan HILIC profiles for a typical mAb. Fluorescence chromatograms for labelled glycans from a mAb using a 2.1 x 150 mm ACQUITY UPLC BEH Amide 130Å 1.7  $\mu$ m column. Peak identifications for the RapiFluor-MS labelled N-glycans are provided. 2-AB labelled N-glycans were prepared using a GlykoPrep® Rapid N-Glycan Preparation with 2-AB kit.

RapiFluor-MS labelled glycans are ideally suited for HILIC chromatographic separations. UHPLC separations with a sub-2- $\mu$ m amide bonded stationary phase such as the 2.1 x 150 mm ACQUITY UPLC BEH Amide 130Å 1.7 $\mu$ m provide the highest sample throughput. **Figure 6** shows UHPLC chromatograms obtained for two samples prepared by different sample preparation techniques, one with 2-AB labelling and the other with RapiFluor-MS labelling.

Notice that each set of labelled N-glycans are resolved by the HILIC separation with comparable selectivities. In this example, the mass load of RapiFluor-MS labelled glycans is lower given the enhanced sensitivity afforded by the label.

Due to the greatly improved MS sensitivity afforded by RapiFluor-MS, HILIC separation of labelled N-glycan samples can now be routinely analysed using LC

combined with fluorescence (FLR) and MS detection. This capability provides high quality data for structural elucidation during characterisation, particularly for low abundance species. In addition, this dramatically improved MS sensitivity enables the opportunity to generate mass data during routine analyses often used in development, production, and quality control environments. Using conventional labelling technologies, the generation of meaningful mass data was only possible with high end mass spectrometers, since traditional reductive amination labels have inadequate ionisation efficiencies to facilitate alternative MS detection. With the introduction of RapiFluor-MS, which yields improved fluorescence sensitivity and dramatically improved MS sensitivity and correspondingly higher charge states in positive ion mode ESI-MS [19], it is now possible to use a simpler, lower cost mass detector for glycan MS analysis. An example of use of this workflow to characterise and routinely interrogate samples is shown in **Figure 7**. In this example, we see that a wide range of N-glycan structures are readily detected with both fluorescence and mass detectors. To further highlight the utility of mass detection, particularly in production environments, separation times can be reduced significantly while incorporating the use of selected ion recording, an intrinsic capability of the mass detector. In this way, critical species can be readily monitored without the need for extensive chromatographic separation. As shown in **Figure 8**, both MAN5 and FA2G1 species can be independently monitored despite the fact that they co-elute in this analysis. The total analysis time is less than 10 minutes, and coupled with rapid labelling, results can be obtained from starting with the intact glycoprotein in less than 40 minutes.

## Conclusion

By combining RapiFluor-MS labelling with rapid deglycosylation and a robust HILIC SPE micro-elution clean-up, an analyst can now complete a reproducible N-glycan sample preparation, from glycoprotein to ready-to-analyse sample, in just 30 minutes. Previous rapid and simple N-glycan sample preparation methods, which took approximately 3 hours to complete, required significant compromises to MS sensitivity. RapiFluor-MS labelling removes this limitation [19], thereby making it possible to routinely use mass detection to expedite accurate mass and MS-MS peak characterisation, or provide greater confidence in routine analyses using a mass detector.

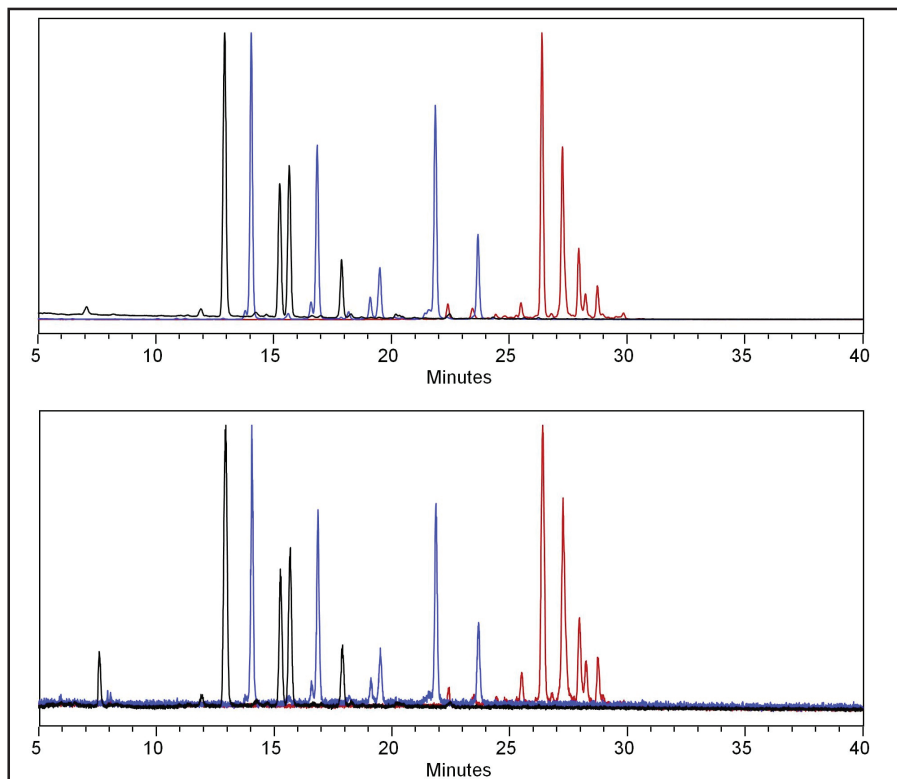


Figure 7. RapiFluor-MS labeled glycans from human IgG (black trace), RNase B (blue trace) and bovine fetuin (red trace) were released with Rapid PNGase F, labelled with RapiFluor-MS reagent. Individual glycan pools were then separated via HILIC and detected with both fluorescence (Top) and mass detection (Bottom).

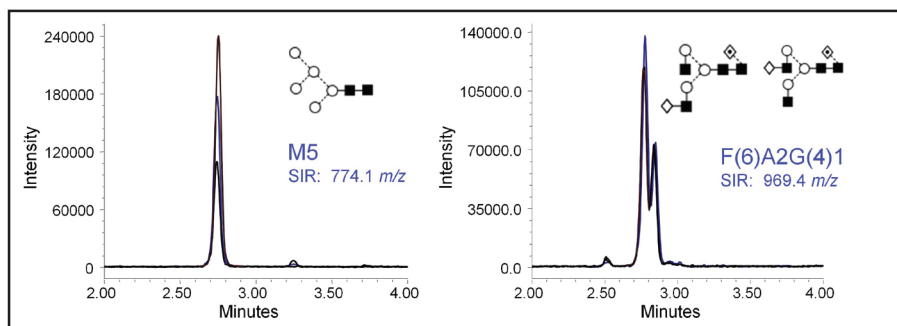


Figure 8. Selected ion recording of RapiFluor-MS labelled MAN5 and FA2G1 species released from IgG separated using HILIC chromatography. Separation was geometrically scaled to a 10 minute analysis time for high throughput.

## References

- Ohtsubo, K.; Marth, J. D., *Cell* **2006**, 126 (5), 855-67.
- Mechref, Y.; Hu, Y.; Garcia, A.; Zhou, S.; Desantos-Garcia, J. L.; Hussein, A., *Bioanalysis* **2012**, 4 (20), 2457-69.
- Ruhaak, L. R.; Miyamoto, S.; Lebrilla, C. B., *Mol Cell Proteomics* **2013**, 12 (4), 846-55.
- Glycosylation Engineering of Biopharmaceuticals*. Humana Press: 2013.
- Rathore, A. S.; Winkle, H., *Nat Biotechnol* **2009**, 27 (1), 26-34.
- Beck, A.; Wagner-Rousset, E.; Ayoub, D.; Van Dorsselaer, A.; Sanglier-Cianferani, S., *Anal Chem* **2013**, 85 (2), 715-36.
- Dalziel, M.; Crispin, M.; Scanlan, C. N.; Zitzmann, N.; Dwek, R. A., *Science* **2014**, 343 (6166), 1235681.
- Tarentino, A. L.; Gomez, C. M.; Plummer, T. H., Jr., *Biochemistry* **1985**, 24 (17), 4665-71.
- Rudd, P. M.; Guile, G. R.; Kuster, B.; Harvey, D. J.; Opdenakker, G.; Dwek, R. A., *Nature* **1997**, 388 (6638), 205-7.
- Tharmalingam, T.; Adamczyk, B.; Doherty, M. A.; Royle, L.; Rudd, P. M., *Glycoconj J* **2013**, 30 (2), 137-46.
- Houel, S.; Hilliard, M.; Yu, Y. Q.; McLoughlin, N.; Martin, S. M.; Rudd, P. M.; Williams, J. P.; Chen, W., *Anal Chem* **2014**, 86 (1), 576-84.
- Kaneshiro, K.; Watanabe, M.; Terasawa, K.; Uchimura, H.; Fukuyama, Y.; Iwamoto, S.; Sato, T. A.; Shimizu, K.; Tsujimoto, G.; Tanaka, K., *Anal Chem* **2012**, 84 (16), 7146-51.
- Ahn, J.; Bones, J.; Yu, Y. Q.; Rudd, P. M.; Gilar, M., *J Chromatogr B Analyt Technol Biomed Life Sci* **2010**, 878 (3-4), 403-8.
- Marino, K.; Bones, J.; Kattla, J. J.; Rudd, P. M., *Nat Chem Biol* **2010**, 6 (10), 713-23.
- Leymarie, N.; Zaia, J., *Anal Chem* **2012**, 84 (7), 3040-8.
- Mechref, Y.; Hu, Y.; Desantos-Garcia, J. L.; Hussein, A.; Tang, H., *Mol Cell Proteomics* **2013**, 12 (4), 874-84.
- Ruhaak, L. R.; Zauner, G.; Huhn, C.; Bruggink, C.; Deelder, A. M.; Wuhrer, M., *Anal Bioanal Chem* **2010**, 397 (8), 3457-81.
- Cook, K. S.; Bullock, K.; Sullivan, T., *Biologicals* **2012**, 40 (2), 109-17.
- Lauber, M. A.; Yu, Y. Q.; Brousniche, D. W.; Hua, Z.; Koza, S. M.; Magnelli, P.; Guthrie, E.; Taron, C. H.; Fountain, K. J., *Anal Chem* **2015**, DOI: 10.1021/acs.analchem.5b00758.